
Installation

Analytical Balances
Analysenwaagen
Balances d'analyse

B5 (H 26)
B6 (H 26)
B6 C 200
B5 D4

METTLER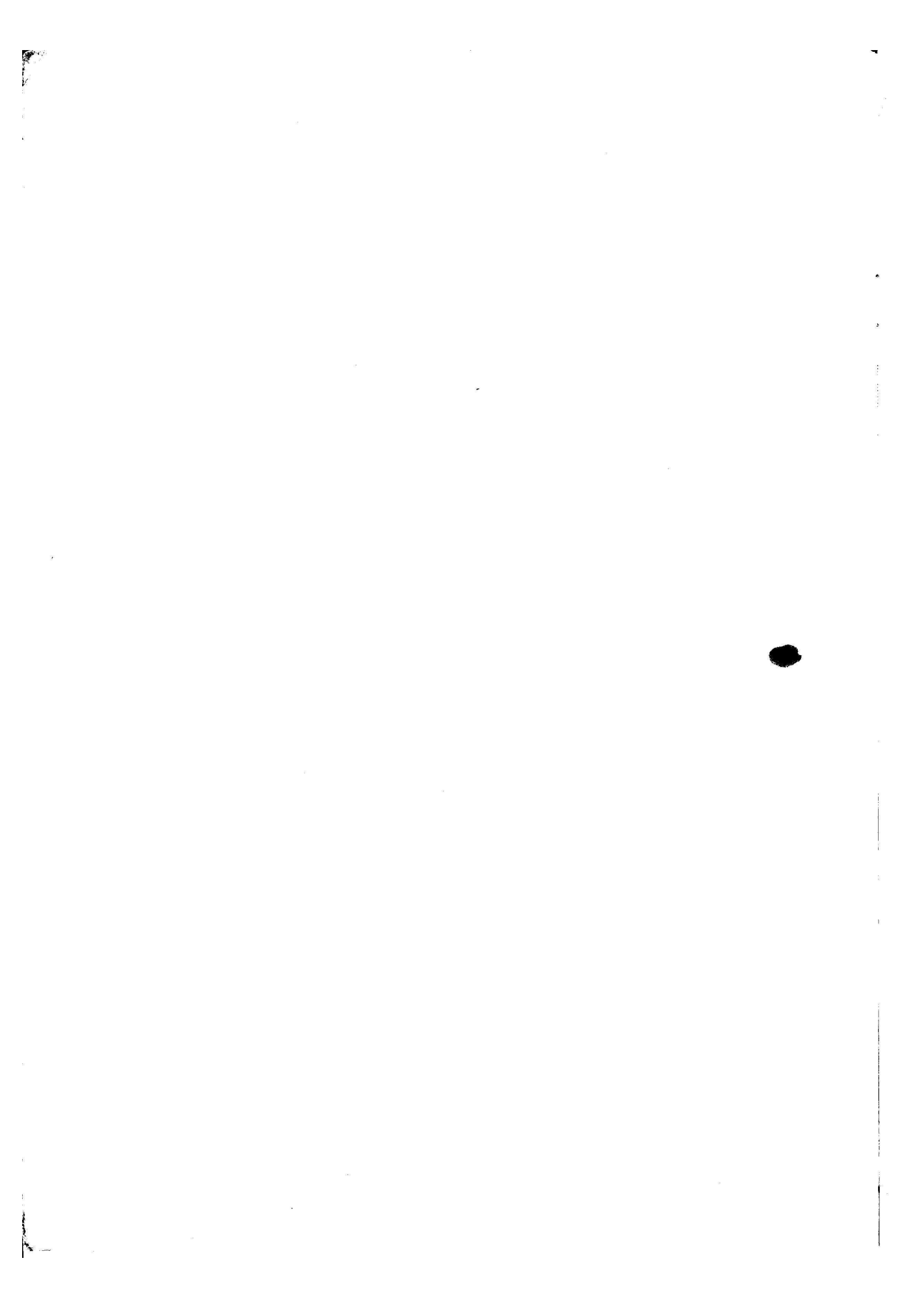

The perfect function of Mettler balances is ensured only if they are set up and operated in accordance with the instructions below.

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Installation Balances B

Contents

I. Unpacking and Setting-up Instructions

- Check voltage
- Scale image sharpness
- Lateral adjustment of scale image

II. Setting Balance to Work

- Zero point adjustment
- Zero point corrections
- Scale deflection (sensitivity) check
- Adjustment of scale deflection

III. Help yourself

- Changing lamp
- Adjusting image brightness
- Minor disturbance

IV. Special equipment

- Taring of suspension
- Weighing below the weighing compartment

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I. Unpacking and Setting-up Instructions

- I. 1. Remove plastic dust cover and packing paper (spares and accessories ordered additionally are enclosed in the wood wool). Remove pads ① and ②.

B...H26:

Remove wooden frame from air freight parcels.

Unpack cable and mains plug.
Place balance where predetermined.

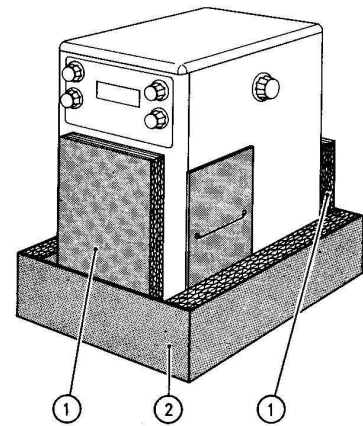
Fig. 1

- I. 2. Remove side members ③ (cut wire). Remove the following parts from the weighing compartment:
 - Plastic cover
 - Two cardboard boxes
 - Weighing pan
 - Bottom glass pad (packing part)
 - Intermediate sheet and protective cardboard

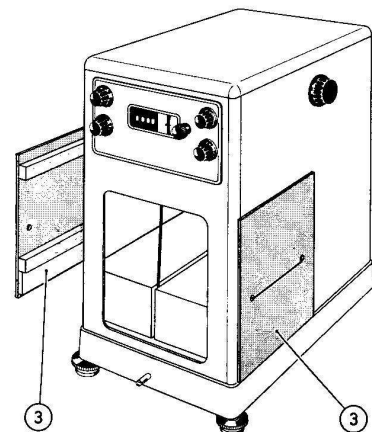
Fig. 2

- I. 3. Lift off housing cover and remove covering board ④.

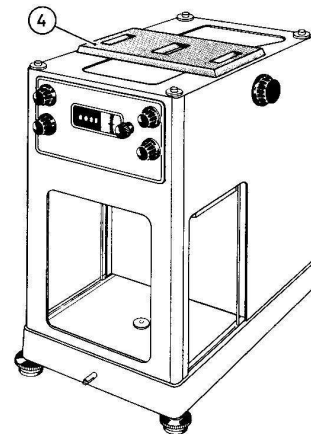
Fig. 3

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- I. 4. Remove weight packing block ⑤ as follows:
 - Loosen both safety nuts ⑥.
 - Turn out both screws ⑦ by a few turns until the built-in weights are released.
 - Hold packing block ⑤ and cut plastic cord.
 - Pull packing block ⑤ downward and remove.

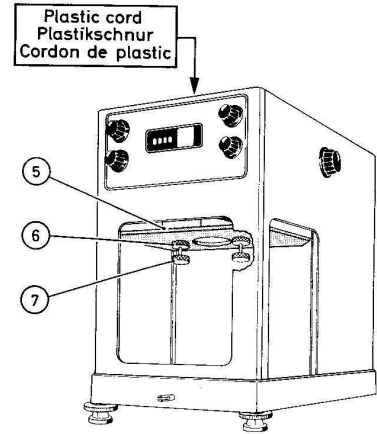
Fig. 4

- I. 5. Unpack tools, transformer plate and accessories from cardboard boxes.

- I. 6. Remove all yellow packing parts in the following order:

- Stirrup ⑭ (four screws).
- Screw ⑫ with rubber disk; place suspension ⑬ carefully forward on the weight levers.
- Cylindrical screw ⑩ and hex head screw ⑧.

Caution: Hold beam and allow it slowly to slide into the bearings.

- Bridge ⑪ (hold bridge) and nylon washer.
- Plate ⑮.
- Both columns ⑨ with washers.

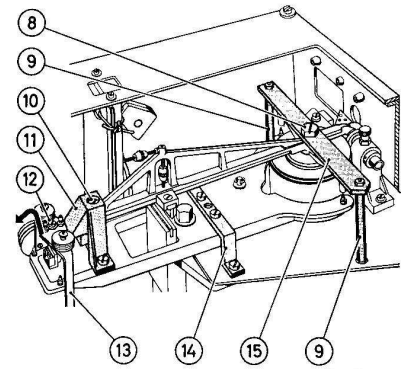
Fig. 5

- I. 7. Remove beam.

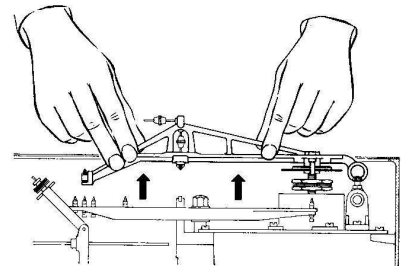
Fig. 6

- I. 8. Check voltage

Remove protection sheets ⑲ and paper pads. Take three screws ⑯ from paper bag. The voltage set on the transformer plate must correspond to that of the supply: if necessary, adjust brass plate ⑳. Attach transformer plate ⑱ to rear panel with the three screws ⑯. Mount protection sheets ⑲.

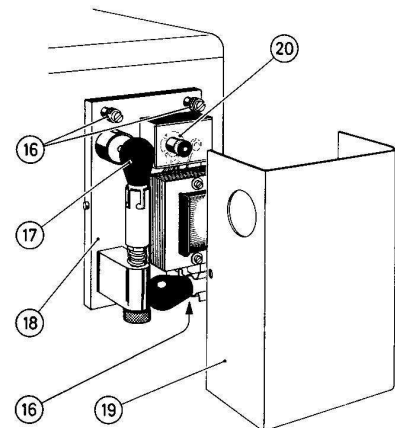
Fig. 7

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Caution: Clean with dry cloth or wash-leather only.

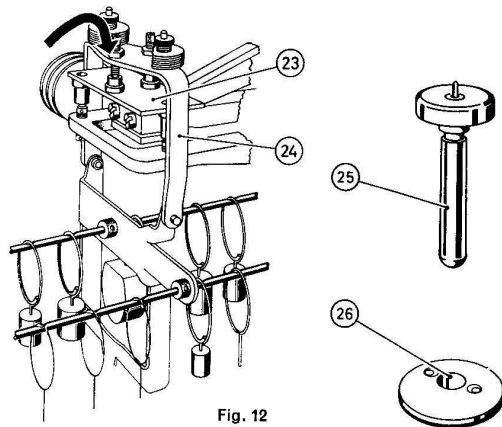
Fig. 12

- I. 12. Insert white intermediate sheet 27 (the two facets facing the front window).

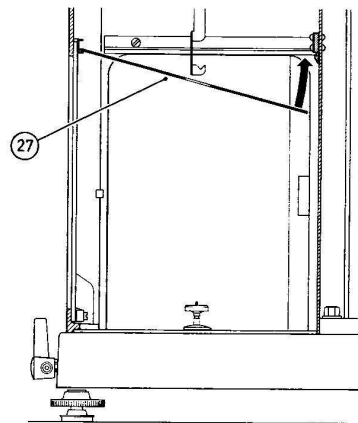
Fig. 13

- I. 13. Clean pan and engage in the pan hook (hook on the pan yoke with opening to the right).

B5 D4:

The pan of these units must be additionally loaded either with the special pan or the compensating weight.

B...GD:

The pan of these units must be additionally loaded either with the hanger or the compensating weight (cf. also IV.2. Weighing below the weighing compartment).

Level balance with the three foot screws. Set weight control to maximum load and slowly return to zero.

Lightly tip pan:

It should come to rest of its own after about four to six oscillations.

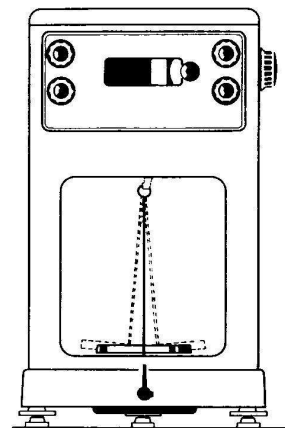
Fig. 14

Effect any necessary corrections as follows:

- Unhook pan.
- Pull pan brake 28 from guide 30.
- The pan brake can be adjusted by turning the sleeve 29.

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Arrest balance (position of arrestment lever, cf. Fig. 20).
 Unload pan and replace housing cover.

Fig. 20

I. 17. Lateral adjustment of scale image

Place housing cover across the balance.

Caution: The two housing clamps must not foul the beam.

Load pan with about 2 g (enclosed hair brush).

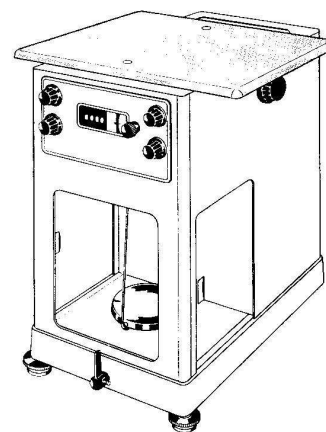
Fig. 21

Semi-release balance (position of arrestment lever, cf. Fig. 22).

Fig. 22

Check overlap (cf. Fig. 24).

Effect necessary corrections as follows:

- Slightly loosen screw ③ and set desired overlap by turning angle prism ②.
- Tighten screw ③ and recheck overlap.

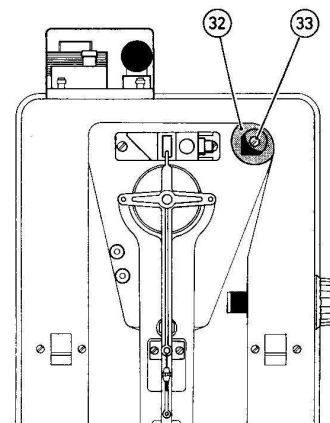
Fig. 23

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B5 D4

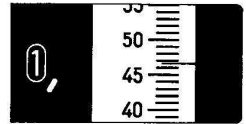
Fig. 24

Arrest balance (position of arrestment lever, cf. Fig. 25).
Unload pan and replace housing cover.

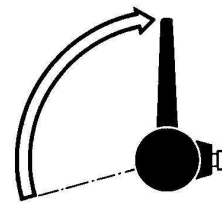
Fig. 25

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II. Setting Balance to Work

Mettler balances are precision instruments of practical and sturdy construction. They are easy to operate if the following rules are observed:

Three fundamental rules for maintaining precision and accuracy

Vibration-free location


It is best to use the balance in a room with constant temperature and humidity. Place the balance on a rigid weighing table which is as heavy as possible (stone table).

Load and unload pan only when balance is arrested

In arrested position, the beam and hanger each rest on three arrestment pins. The knife-edges are fully relieved and will thus not be affected by sudden changes of load such as occur when the pan is loaded and unloaded. The arrestment lever should be moved slowly until the scale lighting is switched on.

The arrestment lever can be shifted on its shaft by pressing on the knob located on the side.

Position A = operating position
Position B = arrestment lever locked

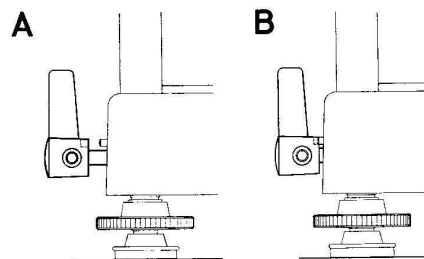Turn weight setting knobs only with balance semi-released

Semi-release protects the knife-edges from damage during weight setting. The sudden load changes resulting from applying and removing weights are to a large extent absorbed by the arrestment pins as the beam movement is limited when semi-released.

Positions of arrestment lever

Balance arrested Position I

Load and unload pan.

Important: Place arrestment lever in position I prior to making any adjustments of the balance.

Balance semi-released Position II

Operate weight control knobs.

Balance fully released Position III

Set zero point.
Read result.

Caution: Check levelling of balance before weighing and adjusting!

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II. 1. Zero point adjustment:

Before checking zero point, make sure that:

- a) Balance is levelled with the three foot screws.
- b) Weight control is set to zero
 - B6; B6 C200; B 6 H26:
Set micrometer to zero.
- c) Side windows are closed.

Fully release balance (arrestment lever in position III).

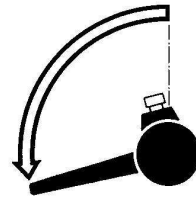
Fig. 26

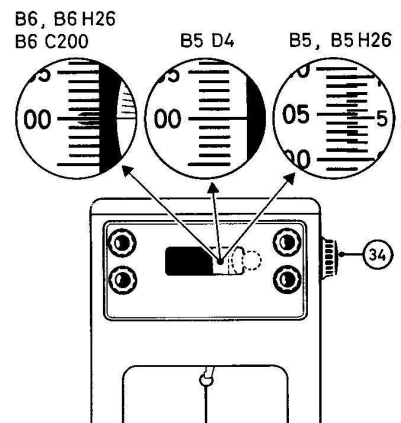
Fig. 27

Turning zero point knob 34 to place zero division of optical scale in the fork and, respectively, at zero line. Effect corrections as per Point II. 2.

Arrest balance (arrestment lever in position I).
Replace housing cover.

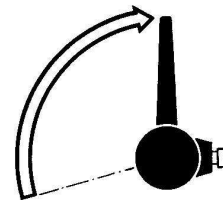
Fig. 28

II. 2. Zero point corrections

If the zero point can be reached with difficulty only or not at all, correction using the fine tare weight 35 will be necessary:

- Set zero adjustment range to center.
- Arrest balance (arrestment lever in position I).
- Remove housing cover.

Correct as per Fig. 29. After correction, check zero point as per Point II. 1.

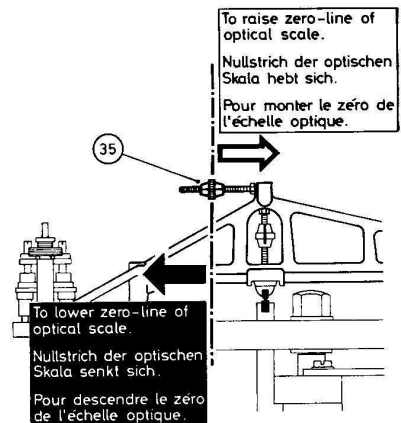
Fig. 29

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II. 3. Sensitivity check

- Set weight control knob to lowest value.
- Load pan with check weight as per Fig. 30.
- Fully release balance (arrestment lever in position III) and set zero point.
- Slowly turn weight control knob back to zero.
- The scale should come to rest exactly at 100. For corrections, if any, see Point II. 4.
- Arrest balance (arrestment lever in position I). Unload pan.

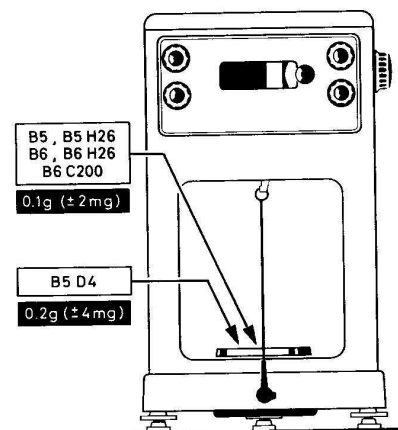
Fig. 30

II. 4. Sensitivity adjustment

If reading is below 100:

- Arrest balance (arrestment lever in position I) and remove housing cover.
- Turn tare weight (36) higher. Replace housing cover and again check sensitivity as per Point II.3.

If reading is above 100:

- Arrest balance (arrestment lever in position I) and remove housing cover.
- Turn tare weight (36) lower, fit housing cover and again check sensitivity as per Point II. 3.

If sensitivity adjustment must be effected using the fine tare weight, repeat sensitivity check (point II. 3.) after about 30 minutes (temperatures balance).

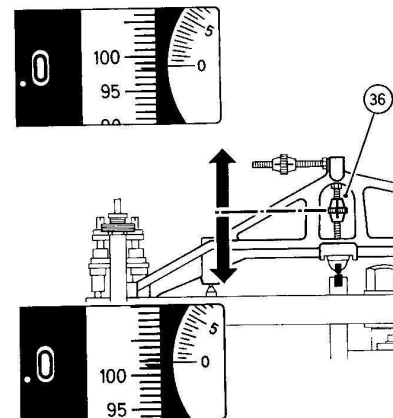
Fig. 31

**The balance is now ready for operation.
For weighing instructions, see separate
weighing sheet.**

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III. Help yourself

Changing lamp

Remove protection sheet and replace defective or dim lamp.
Check image brightness.

Important: Order another spare lamp immediately!

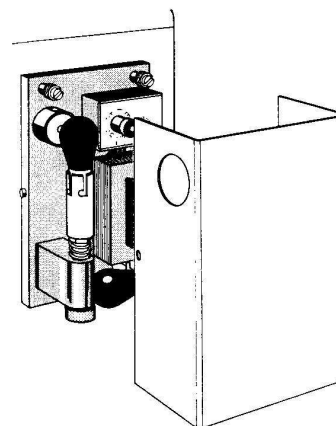
Fig. 32

Adjusting image brightness

Semi-release balance.
Adjust image brightness by vertical adjustment of the lamp.

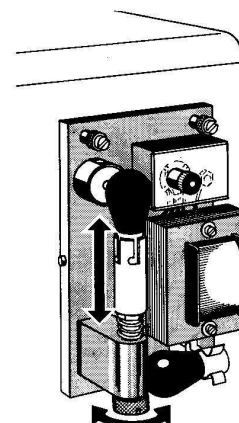
Fig. 33

Minor disturbance

If the scale moves jerkily or blocks after releasing:

- Arrest balance.
- Turn weight control knobs to maximum load and slowly back to zero.
- Clean pan brake (37) and pan brake guide (38).

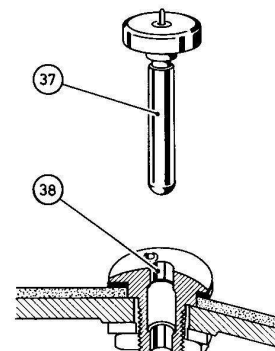
Fig. 34

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IV. Special equipment

IV. 1. Taring of suspension

If a weight difference in the suspension must be compensated, e. g. after installing pan hooks, this is effected by removing taring disks [41].

- Weight of large disks: about 1 g
- Weight of small disk: about 0.1 g

Important: Remove the same number of disks on both the left and the right.

Fine corrections are performed by means of the fine tare weight (40).
Removed taring disks should be placed on the screws (39).

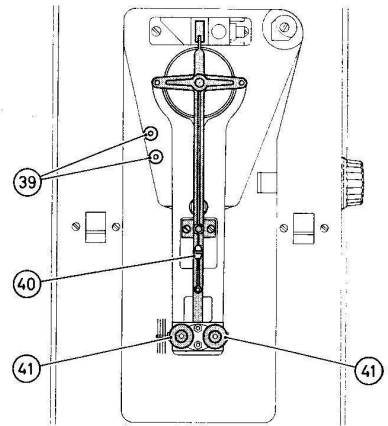
Fig. 35

IV. 2. Weighing below the weighing compartment

For weighings **below** the weighing compartment, the pan must be provided with the hanger.

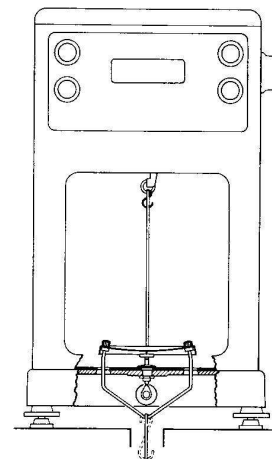
Fig. 36

For weighings **in** the weighing compartment, the hanger is removed and replaced by the extra weight. The weight of the latter corresponds to that of the hanger.

For weighings **in** the weighing chamber, the GD openings can be closed by the enclosed stops.

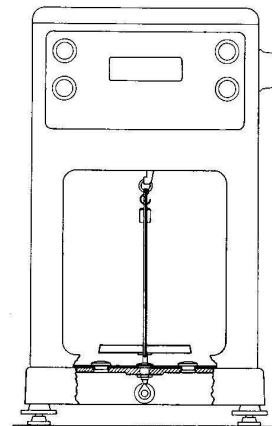
Fig. 37

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